

Bench-Top Capsule Filling Systems

For Use in R&D • Pharmacy Compounding • Nutritional Supplements • Clinical Studies





ProFiller 3700 is the new standard in 300-hole bench-top capsule filling systems. Available with a European style Semi-automatic Orienter as ProFiller 3800. Experienced users like the higher output, fast size changes, easy to clean design & optional accessories. First time users love the ease-of-use & upgrade options.

Easier To Use

- Easier to use than other 200 or 300 hole systems.
- Indicator gauges to assist in capsule orientation, separation and locking.
- Shipped fully assembled. Ready for use in less than 1 hour.
- Video tips at <u>www.profiller.com</u>. Well-illustrated manual in multiple languages.

Higher Output

- Unique features include lightweight, 2-piece Orienter (3700), Orienter locater and oversize Powder Tray & Push Plate.
- Tamper size range optimized for higher fill weights and faster filling.
- Adjustable capsule body height to make over-filling easier & faster.
- Fully locks capsules after filling.

Pharmacopoeia Compliant

- Fill weights, USP/EP method within ± 1-3% of average. Typically all 300 capsules/batch are within ± 4-7% of target. More information at www.profiller.com/3fw.
- Validation package available for cGMP compliance.
- Works with all capsules including gelatin, vegetarian/HPMC and starch capsules.

The easiest 300-hole filler to use! Up to 6000 to 9000* capsules/hour.

*based on powder, operator speed & technique, excipients and weight variation limits. Output is similar for ProFiller 3700 and 3800.

Grows With You

- Add a Vibrator at any time.
- Change Parts for 3700 in Size 000-5, 00el-2el & AAA-E; for 3800 in Size 00-4.
- Order Filler without a Tamper for vibration only operation or in a reduced space configuration (RS) for use under smaller powder hoods.

GMP & Faster Clean-Out

- Dishwasher safe Inox (SS316) powder contact parts.
- Designed for fast disassembly & visual verification of cleaning.
- Filler base in lightweight, corrosion resistant anodized aluminum or upgrade to optional Inox (SS316).
- Filler's gear, cam and optional Vibrator are fully covered.

Exclusive Accessories


- Partial Batch Kit for filling fewer than 300 capsules per batch.
- Capsule Counter.
- Production Stand.
- Over-Encapsulation attachments for tablets, capsules & caplets
- Caps Tray in Inox (SS316) or optional lightweight anodized aluminum.

Distributed by:

CAPSUGEL

Toll Free Phone: 888-783-6361

www.capsugel.com

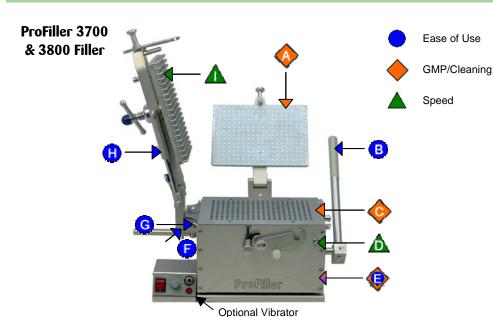



SS316 (Inox) Locking Plate. No rubber pad to wear out or clean



Rear locking handle for fast, easy movement across front of machine. New internal gear design for easier locking



Dishwasher safe Inox (SS316) powder contact parts including Bottom Plate (not visible)



Adjustable capsule body height for easier & faster over-filling of capsules



Base Frame in light-weight anodized aluminum or upgrade to Inox (SS316)



Cam position indicator, great for training new operators



Over-size thumb screws for easy size change over




Dual springs make Tamper easier to lower



Tamper pin size range optimized for faster filling & higher fill weights. Size ranges: 000 / 00 / 0-1 / 2-3 / 4 / 5



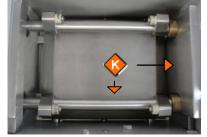







Cover for gear mechanism & Vibrator to minimize powder entry (inside view)




Powder Tray is double-depth for faster Spreading




Multi-use Push Plate speeds up operations

3700 Orienter

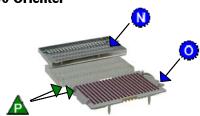

3800 Orienter

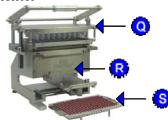

- N
- 2-piece Hand-Held Orienter. Shake only the lightweight Orienter Tray
- 0
- capsules are fully locked Locating holes in Orienter and tapered locating pins for fast & accurate

orientation of capsules into Caps Tray

Locking indicator instantly confirms

- Q
 - Orienter Base common for Size 00-4. Traditional European style Orienter for operators who prefer it
- Guide lines to locate Caps Tray for partial R batches
 - Locking indicator instantly confirms capsules are fully locked

Adjustment-Free Size Changeovers

Shipped Fully Assembled





ProFiller 3700 Ordering Information

For Use in R&D • Pharmacy Compounding • Nutritional Supplements • Clinical Studies

ProFiller 3700 Bench-Top Filler & Hand-Held Orienter



ProFiller 3700		
Capsule Size	Item Number	
00	P7T BHS 00	
0	P7T BHS 0	
1	P7T BHS 1	
2	P7T BHS 2	
3	P7T BHS 3	
4	P7T BHS 4	
5	P7T BHS 5	
000*	P7T BHS 000	
For 00el*, 0el	- 2el, AAA-E sizes contact us.	

^{*}Standard 180 hole Filler (192 holes special order).

ProFiller 3700 System includes everything you need to fill one capsule size. The standard configuration is shown above with the optional Vibrator. The standard Bench-Top filler has a Tamper and a lightweight anodized aluminum Base Frame. For alternate configurations including an Inox (SS316) Base Frame see the FAQ page.

Included at no extra charge: a Video CD, illustrated manual, plastic storage box for easy storage of loose items when the machine is not in use, empty and locked Capsule Bins, and a set of empty capsule Measuring Bowls.



ProFiller Bench-top Vibrator


Optional Vibrator recommended for faster filling of powders that bridge or are granular, fluid or flour-like. It also reduces weight variation of fine powders. Fully enclosed in Inox (SS316). Filler sits over Vibrator. Also available with Inox (SS316) Base Plate.



Item #	Voltage/ Frequency
P78 VIB 110	110-120V/50Hz
P78 VIB 230	220-240/50 Hz

ProFiller 3700 Change Parts

	Change Parts	Tamper	
Capsule Size	Item #	Item #	
00*	P7T CPS 00	P78 T300P 00	
0*	P7T CPS 0	P78 T300P 0-1	
1*	P7T CPS 1	P78 T300P 0-1	
2*	P7T CPS 2	P78 T300P 2-3	
3*	P7T CPS 3	P78 T300P 2-3	
4**	P7T CPS 4	P78 T300P 4	
000, 5, 00el - 2el & AAA-E	visit www.profiller.com/3700cps		

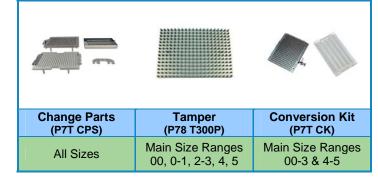

^{*} If you have a 3700 system in Size 4 only, also order Conversion Kit P7T CK 00-3.

You also need to order Conversion Kit P7T CK 4-5, unless you can already fill Size 5 on your system. All Change Parts include a plastic storage box that can also hold a Tamper. Other sizes available upon request.






ProFiller 3800 Ordering Information

For Use in R&D • Pharmacy Compounding • Nutritional Supplements • Clinical Studies

ProFiller 3800 Bench-Top Filler & European Style Orienter

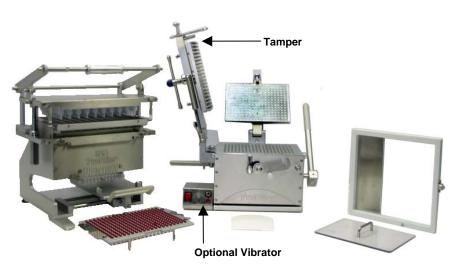

ProFiller 3800			
Capsule Size	Item Number		
00	P8S BSA 00		
0	P8S BSA 0		
1	P8S BSA 1		
2	P8S BSA 2		
3	P8S BSA 3		
4	P8S BSA 4		
Sizes 000, 5, 00el - 2el, AAA-E available in ProFiller 3700.			

ProFiller 3800 System includes everything you need to fill one capsule size. The standard configuration is shown above with the optional Vibrator. The standard Bench-Top filler has a Tamper and a lightweight anodized aluminum Base Frame. For alternate configurations including an Inox (SS316) Base Frame see the FAQ page.

Included at no extra charge: a Video CD, illustrated manual, plastic storage box for easy storage of loose items when the machine is not in use, empty and locked Capsule Bins, and a set of empty capsule Measuring Bowls.



ProFiller Bench-top Vibrator


Optional Vibrator recommended for faster filling of powders that bridge or are granular, fluid or flour-like. It also reduces weight variation of fine powders. Fully enclosed in Inox (SS316). Filler sits over vibrator. Also available with Inox (SS316) Base Plate



Item #	Voltage/ Frequency		
P78 VIB 110	110-120V/50Hz		
P78 VIB 230	220-240/50 Hz		

ProFiller 3800 Change Parts

	Change Parts	Tamper	
Capsule Size	Item #	Item #	
00*	P8S CPS 00	P78 T300P 00	
0*	P8S CPS 0	P78 T300P 0-1	
1*	P8S CPS 1	P78 T300P 0-1	
2*	P8S CPS 2	P78 T300P 2-3	
3*	P8S CPS 3	P78 T300P 2-3	
4**	P8S CPS 4	P78 T300P 4	
000, 5, 00el - 2el & AAA-E	visit <u>www.profiller.com/3700cps</u>		

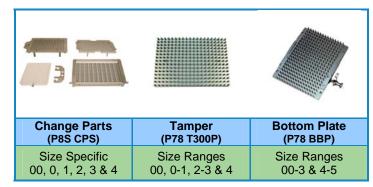

^{*} If you have a 3800 system in Size 4 only, also order Bottom Plate P78 BBP 00-3.

^{**} You also need to order Bottom Plate P78 BBP 4-5, unless you can already fill Size 5 on your system using a hand-held orienter. All Change Parts include a plastic storage box.




ProFiller Inox (SS316) Ordering Information

For Use in R&D • Pharmacy Compounding • Nutritional Supplements • Clinical Studies

ProFiller Stainless Steel Upgrade

For high GMP requirements the ProFiller 3700 and 3800 can be ordered in a Stainless Steel 316 (Inox) version. All Materials of Construction are the same as the standard ProFiller except for those shown here which are in Inox (SS316).

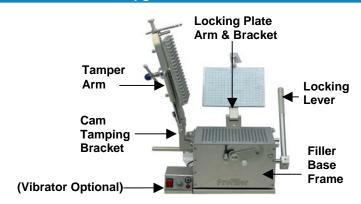



ProFiller 3700 Inox (SS316)

Capsule Size	Item Number	
00	P7T BHS 00 SS	
0	P7T BHS 0 SS	
1	P7T BHS 1 SS	
2	P7T BHS 2 SS	
3	P7T BHS 3 SS	
4	P7T BHS 4 SS	
5	P7T BHS 5 SS	
000*	P7T BHS 6 SS	
00el-2el, AAA-E sizes contact us		



ProFiller 3800 Inox (SS316)

Capsule Size	Item Number	
00	P8S BSA 00 SS	
0	P8S BSA 0 SS	
1	P8S BSA 1 SS	
2	P8S BSA 2 SS	
3	P8S BSA 3 SS	
4	P8S BSA 4 SS	
000, 5, 00el*, 0el - 2el, AAA-E only available for 3700.		

Stainless Steel Vibrator

Heavy duty Vibrator for use ProFiller 3700/3800 in Inox (SS316). Includes full Inox (SS316) base.



Item #	Voltage/ Frequency
P78 VIB 110 SS	110-120V/50Hz
P78 VIB 230 SS	220-240/50 Hz

Change Parts

Change Parts are all Inox (SS316).

To order Change Parts for ProFiller 3700 or 3800 lnox (SS316) see page 3 or 4.



^{*}Sizes 000 and 00el in 180-holes (192 holes special order)



Bench-Top Capsule Filling System

For Use in R&D • Pharmacy Compounding • Nutritional Supplements • Clinical Studies

Installation Requirements




Width:


48 inches (120cm) is recommended width if Filler and Orienter are operated on same table.
34 inches (90 cm) is minimum width to operate Filler if Orienter is kept on another table.

Depth: 30 inches (75cm) depth is recommended. 24 inches (65cm) is the minimum depth.

Height: 29 inches (73cm) or office desk height is recommended.


Electricity is required for optional Vibrator. No compressed air or water is required for operation.

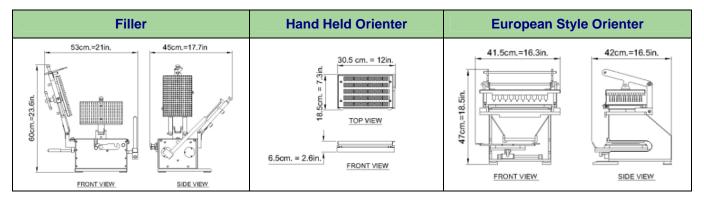

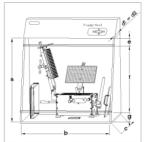

For detailed measurements and information on selecting a powder hood, visit http://www.profiller.com/powder-hood.

Shipping Information

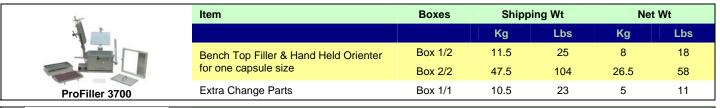





For Inox (SS316) Filler on page 5, add 9.5 kg (28 lbs) to shipping and net weight of Box 2. If optional Inox (SS316) Vibrator ordered, add an additional 13 kg (29 lb) to shipping and net weight.

Inox (SS316) Filler

For ProFiller 3700 & 3800 with Aluminum Vibrator add 10 kg (22 lb) weight to shipping weight as well as net weight of Box 2.





Optional Accessories

For Use in R&D • Pharmacy Compounding • Nutritional Supplements • Clinical Studies

Opti	onal Accessories (ProFille	r 3700 / 3800)		
Item Description		Voltage/Frequency	Capsule Size*	Item #
Powder Vibrator *		110-120 Volts/60Hz	All Sizes	P78 VIB 110
Recommended for faster filling of powders that bridge or are granular, fluid or flour-like. Also reduces weight variation of fine powders. Fully enclosed in lnox (SS316). Filler sits over Vibrator. Powder Tray Clamps included.	9	220-240 Volts/50Hz	All Sizes	P78 VIB 220
Powder Tray Clamps For pharmacy compounding and very fine powders. Reduce powder leakage under the Powder Tray. Available October 2011	21	7	All Sizes	TBD
Production Stand Reduce the space required to operate the ProFiller. The stand organizes and provides easy access to all the loose parts. Stand only. (Parts shown for illustration)			All Sizes	PXA STAND
Lifting Handle Attachment Attaches to Lifting Handle to make capsule separation and locking under a powder hood easier.			All Sizes	P78 LH PH
Over Fill (OF) Change Parts Standard system allows capsule bodies to be lowered 2 mm. OF Change Parts allow for an additional 2mm of space for twice the overfill capacity.			000-1	TBD †
Capsule Counters Divide your batch of capsules into 10 groups of 30 or 3 groups of 100 capsules.	90 e.e.	Anodized Aluminum	All Sizes	PXA CTR 3X100 PXA CTR 30X10
Opti	onal Accessories (ProFiller	r 3700 ONLY)		
Capsule Locking Plate & Capsule Locker For fully locking capsules with coarse or sticky powders e.g. salt, dolomite, frankincense, charcoal.	17		00-4	PXA CLP 00-4
Capsule Locking Plate is common for sizes 00-4. Both Locker & Locking Plate are required. Capsule Locker is size specific and locks 25 capsules at a time.			00-4	PXA CL [†]
Partial Batch Kit Fill fewer than 300 capsules per batch. Includes 1 each of Orienter Divider, Powder Tray Divider & Mini Powder Spreader. Order 30-pin Tamper separately.	SPECIAL SEC		All Sizes	P67 PBK
30-pin Tamper For use with ProFiller 3700 Partial Batch Kit. Tamper in Sizes 00, 0-1, 2-3, 4.			00-4	P67 T30P †
Caps Tray For 3700 Purchase an additional Caps Tray		Anodized Aluminum**	All Sizes	P67 CT AL †
**Light-weight Anodized Aluminum Caps Tray is not Dishwasher Safe.		Inox (316 SS)	All Sizes	P67 CT SS †

[†] Specify Capsule Size when ordering, some accessories are not available for special sizes (000, 5, 00el-2el & AAA-E). * ProFiller 3700 & 3800 can be ordered without the Tamper Assembly for use only on the Vibrator. See page 8.



Optional Accessories

For Use in R&D • Pharmacy Compounding • Nutritional Supplements • Clinical Studies

Optional Accessories (ProFiller 3800 ONLY)				
Item Description			Capsule Size	Item #
Caps Tray For 3800	anninininininininininininininininininin	Anodized	00-4	P8 CT AL [†]
Purchase an additional Caps Tray		Aluminum*		
*Light-weight Anodized Aluminum Caps Tray is not Dishwasher Safe.		Inox (316 SS)	00-4	P8 CT SS [†]

Over-Encapsulation Accessories (ProFiller 3700 / 3800)					
Item Description		Capsule Size	Item #		
Tablet Fill Attachment Over-encapsulate round tablets 6.0-7.3 mm diameter	A	Standard	PXA TFA STD		
and 2.0-3.5 mm height. For round tablets outside this size range, order a custom TFA (delivery 4-8 weeks).		Custom	Tablet Dimensions required		
Orienter for Capsule-in-Capsule		00-4	P67 NEC OCPOB [†]		
This allows you to put capsules inside capsules using your existing ProFiller System. Order OE Locator Tray separately.		0el – 2el	P67 NEC OCPOB [†]		
Caplet Attachment					
For caplets or non-round tablets. Engineering drawing of tablet required. Delivery 4-8 weeks. Order OE Locator Tray separately.		00-4, 0el – 2el	Caplet Dimensions Required		
	Control of the Contro				
OE Locator Tray		All Sizes	PXA OFF		
Required for use with Orienter for Capsule in Capsule or Caplet Attachment.		All Sizes	PAA OEF		

 $^{^{\}dagger}$ Specify capsule size

3700/3800 Filler Base Options









	High GMP Model ^a	No Tamper ^b	Configurations For Low Height Powder Hoods Video: www.profiller.com/llpvideo	
	Base Frame in Inox (SS316) See p. 5	Filler without Tamper. Sold only with Vibrator.	Filler with loose Locking Plate & special Handle. Tamper is on rear of Filler.	Filler with loose Locking Plate, special Handle & no Tamper. Sold only with Vibrator.
Item Code Suffix		add NT	add SLP	add SLPNT

 $^{^{\}rm a}$ High GMP model is also available in any of the configurations shown above. $^{\rm b}$ Tamper arm & Tamper (in any one size) can be ordered later for \$950.





Frequently Asked Questions

For Use in R&D • Pharmacy Compounding • Nutritional Supplements • Clinical Studies

Frequently Asked Questions

1. What is the output?

ProFiller 3700 and 3800 fill up to 6000 – 9000 capsules per hour for most powders. Our engineers have included features, some patented, that make the ProFillers faster and easier to use. The lower output is for fine powders. The higher output is for coarse free flowing powders. Your results will vary based on powder flow properties, operator speed & technique and use of excipients.

2. What types of capsules can I fill?

ProFillers have been tested with different brands and types of capsules including gelatin, vegetarian/HPMC and starch capsules.

3. What size capsules can I fill?

The 3700 fills sizes 000-5, 00el-2el, and AAA-E. The 3800 fills sizes 00-4. To order Change Parts for the 3700 see page 3 & for 3800 see page 4.

4. Do you have more information on Fill Weight Variation & Achieving Higher Fill Weights?

Fill weights, USP/EP method, within \pm 1-3% of average. When weighed individually (a stricter standard than USP/EP) typically all 300 capsules/batch are within \pm 4-7% of target. For more information visit www.profiller.com/3fw.

5. Why Might I Need a Conversion Kit?

The Filler's Bottom Plate (3700/3800) and the Hand-held Orienter Base (3700) come in different capsule size ranges: 000, 00/0/1/2/3 and 4/5. For capsule sizes outside the size range of your system, you need a Conversion Kit along with Change Parts.

6. Can I see the system before I buy?

You can view our online videos at www.profiller.com

7. Can I over-encapsulate or double-blind tablets, capsules and caplets?

Yes, you can. See page 8 and then contact us for more information.

8. Can I fill liquids or sticky powders?

Yes, you can fill capsules with sticky powders or non-aqueous liquids. Capsules filled with liquids for stability or clinical trials, if not sealed, will need to be kept vertical to minimize leakage between the cap and body. Capsules filled with non-aqueous liquids for commercial sale require sealing using equipment purchased separately. When filling sticky powders, use of a Vibrator can improve performance by eliminating Tamper pin contact with powder. For more information visit www.profiller.com/vibrators/use.

9. Can the ProFiller be cleaned in an automatic dishwasher or autoclaved?

All powder contact parts are dishwasher safe to 70 °C (160 °F). The SS316 (Inox) Base Frame is autoclave safe. For more information visit www.profiller.com and download the manual.

10. What are the Materials of Construction?

For more information visit www.profiller.com and download the manual.

11. What are the installation and utility requirements?

See page 6 for the recommended table size and height and information on operating under a powder hood. No utilities (electricity, compressed air or water) are required unless you order the optional Vibrator. The Vibrator operates on 110-120 Volts / 60 Hz or 220-240 Volts / 50 Hz. If you have different electrical requirements specify them in your order.



ProFiller Products

For Use in R&D • Pharmacy Compounding • Nutritional Supplements • Clinical Studies

Equipment for Sizes 000 - 5



Handheld 100-hole Capsule Filler with Orienter

Up to 2000-3000 Capsules per hour. Easy to operate and up to 35% faster than similar fillers due to its patented design. Change Parts and a full range of accessories are available. Offered in Sizes 000-5 and all el sizes. Over 5000 in use globally! ProFill OE for DBcaps[®].





Handheld 300-hole Capsule Filler

For the growing company and for use under powder hoods or in glove-box isolators. Fully locks capsules after filling. The only 300-hole hand-held you can upgrade as the Change Parts of the 3600 also fit the 3700 bench-top Filler. The same Orienter works with 3600 & 3700. Output up to 4500-6900 Capsules per hour. Available in Sizes 000-4 & all el sizes.





Bench-top* Capsule Filler & Hand-held Orienter

Easiest bench-top system to set-up and use. Includes a hand-held Orienter. Up to 6000-9000 capsules per hour. Options include Partial Batch Kit and Vibrator. Available in Sizes 000-5, all el sizes & for DBcaps[®].





Bench-top* Capsule Filler & European Style Orienter

Our bench-top system for those who prefer the traditional European style semi-automatic Orienter. Up to 6000-9000 capsules per hour. Optional Vibrator. Available in Sizes 00-4. Other sizes use a handheld orienter.



*Bench-top Systems are designed for easier operation to enable continuous production. The 3700 includes our innovative hand-held orienter. For customers who use the traditional European style "semi-automatic" orienter we offer the 3800.

Other Equipment



Capsule / Caplet Film Coater

For R&D. Ideal for enteric or film coating small batches of capsules and caplets with solvent based coatings.





Filling Funnels

For precision capsule filling on a balance. Includes Tamper. Made from Inox (316 stainless steel). Available in all capsule sizes including veterinary sizes.



Veterinary Capsule Fillers



ProFill Vet

Handheld filler for sizes 13-Su07. Up to 800 capsules per hour. Lightweight design for laboratory and test-market trials.



Veterinary Capsule System

Bench-top filler for sizes 13-Su07. Up to 1500-2400 capsules per hour with Orienter. Heavy-duty design for production use.

